



ExoGlow-NTA™ Fluorescent Labeling Kit (for Malvern NanoSight)

Cat# EXONTA200A-1

User Manual

See List of Components for storage conditions

Version 3
9/23/2022

A limited-use label license covers this product. By use of this product, you accept the terms and conditions outlined in the License and Warranty Statement contained in this user manual.

Contents

Product Description.....	2
List of Components.....	2
Storage.....	2
General Information.....	2
Protocol for sample labeling.....	2
Protocol to remove free dye	3
Example Data and Applications.....	4
Technical Support.....	5
Licensing and Warranty Statement	6

Product Description

SBI's ExoGlow-NTA™ Fluorescent Labeling Kit (for Malvern NanoSight) (Cat #EXONTA200A-1) enables highly specific quantitation of extracellular vesicles (EVs) from a wide variety of biofluids and isolation protocols. Compatible with Malvern's NanoSight range of NanoTracking Analysis (NTA) instruments (equipped with 488nm laser) to measure size and concentration of EVs, the ExoGlow-NTA kit leverages the fluorescent capabilities of the platform to specifically detect only labeled EVs in heterogeneous mixtures. The proprietary dye formulation in our ExoGlow-NTA kit binds specifically and efficiently to the membranes of intact EVs only, thus avoiding detection of protein aggregates, membrane fragments and other background particles and ensuring much more accurate EV particle counts than currently available methods.

The kits come with four components: 1) Labeling dye 2) Standards 3) Reaction Buffer, and 4) Size exclusion columns. Simply mix the dye with the reaction buffer, add 1-100ug of EVs (or protein equivalent), incubate for 3-5 minutes, and you are ready for fluorescent NTA analysis. The provided standard is a size-controlled synthetic liposome prep that provide a positive control for the NTA particle tracking as well as labeling using the ExoGlow-NTA kit.

List of Components

Item	Volume	Storage Temperature
Reaction Buffer	160 µl	Room Temperature
Labeling Dye	60 µl	Room Temperature
Standards	3 µl	4°C
Size exclusion column	13	Room Temperature

*The kit is for 10 labeling reactions. We will provide 3 reactions for the Standard, as it will not be necessary to run the standard every time the NTA is run.

Storage

The kit is shipped on ice and should be **stored** as indicated in the table above. Properly stored kits are stable for 6 months from the date received.

General Information

The reaction size is based on using 1-100 µg of total proteins in the sample.

We recommend to pre-warm Reaction Buffer for 3-5 min at 37°C and mix well before use. Protect labeling dye from light.

Protocol for sample labeling:

1. Pre-warm Reaction buffer for 3-5 minutes at 37°C and mix well.
2. For each Sample and Standard, add 4 µl of Labeling Dye into 12 µl of Reaction Buffer and mix well until dye is dissolved completely to make Labeling Reaction buffers.

3. Add exosomes (equivalent to 1-100 µg of protein) to the Sample from step 2 and bring total volume of the reaction up to 50 µl with water or 1xPBS. To label Standards, add 1 µl of the standards into Labeling Reaction buffer.

Compound	Sample Reaction	Standard Reaction
Reaction Buffer	12 µl	12 µl
Labeling dye	4 µl	4 µl
Exosomes (1-100 µg)	X µl	-
Standards	-	1 µl
Water or PBSx1 (filtered)	Up to 50 µl	Up to 50 µl

3. Mix well by pipetting and incubate for 3-5 minutes at RT. The tubes do not need to be rotated during the incubation period.

! Protect the tubes from light.

Protocol to remove free dye:

1. Mix the medium inside the column by vortexing.
2. Loosen the cap and twist off the bottom closure.
3. Place column into a collection tube and remove the storage solution by centrifugation at 800 x g for 1 min.
4. To wash the column, add 400 µL filtered PBS and centrifuge at 800 x g for 1 min.
5. Discard flow-through and replace the collection tube.
6. Repeats steps 4 and 5 four more times.
7. Replace the used collection tube with a new clean 1.5 ml eppendorf tube (not provided) for sample collection.
8. Adjust the volume of the sample to 140 µL with 1x1XPBS. Apply 140-180 µL of your sample slowly in the middle of the packed bed of the column. Recommended sample volume is 140 µL.
9. Centrifuge at 800 x g for 1 minute to collect labeled EVs.
10. The samples are ready for fluorescent NTA analysis.

We recommend optimizing your instrument settings with labeled standards for fNTA analysis. The following setting parameters are provided as suggested starting points for instrument optimization:

Instrument	Camera	Camera level	Slide shutter	Slide gain	Shutter/ms	Frames rate/fps	Syringe pump speed
LM 10/ NS 300	CMOS	11-16	890- 1,300	146-512	22.25-32.5	24.98	100

Example Data and Applications

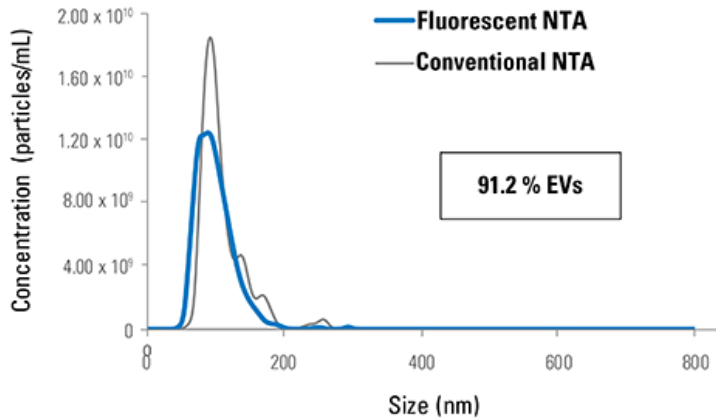


Figure 1. ExoGlow-NTA-labeled liposomes deliver consistent NTA data whether in light scattering or fluorescent mode. The high concordance of NTA and fluorescent NTA data collected from the ExoGlow-NTA Kit internal standards (ExoGlow-NTA-labeled synthetic liposomes) demonstrates the labeling efficiency of the ExoGlow-NTA Dye and accuracy of the fluorescent NTA method for characterizing EVs. **Data collected using NanoSight LM10 instrument equipped with 488nm laser.**



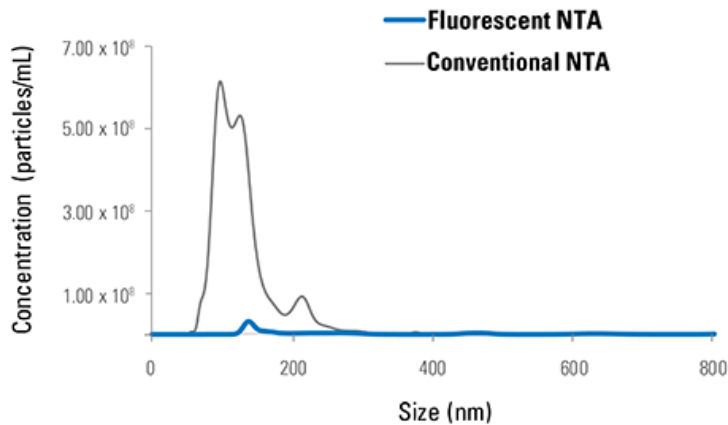


Figure 2. ExoGlow-NTA delivers undetectable background signal. When analyzing the ExoGlow-NTA dye alone in PBS, conventional NTA picks up background particles in the absence of EVs, while fluorescent NTA of the ExoGlow-NTA dye alone shows bias-free undetectable autofluorescence. **Data collected using NanoSight LM10 instrument equipped with 488nm laser.**

A. ExoQuick

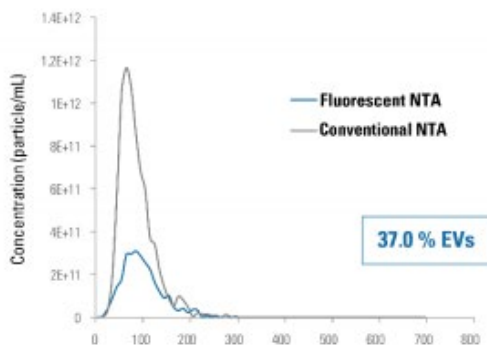

B. Ultracentrifugation + Wash

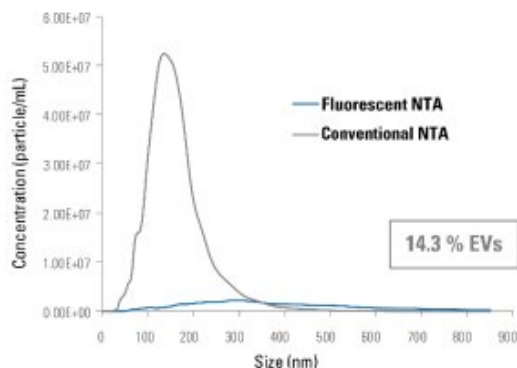

C. Column-based

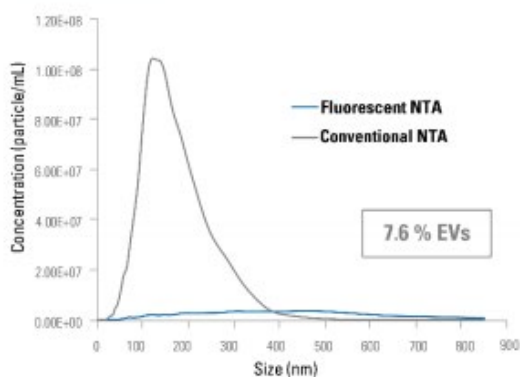


Figure 3. ExoGlow-NTA demonstrates that conventional NTA overestimates EV concentration in samples irrespective of EV isolation method. Representative data comparing conventional NTA and fluorescent NTA for EVs isolated using (A) ExoQuick, (B) ultracentrifugation and wash, or (C) column-based isolation, shows just how much of the conventional NTA signal is due to non-EV particles. ***Data collected using NanoSight LM10 instrument equipped with 488nm laser.***

Technical Support

For more information about SBI products and to download manuals in PDF format, please visit our web site: <http://www.systembio.com>

For additional information or technical assistance, please call or email us at:

System Biosciences (SBI)
2438 Embarcadero Way
Palo Alto, CA 94303

Phone: (650) 968-2200
Toll-Free: (888) 266-5066
Fax: (650) 968-2277

E-mail:

General Information: info@systembio.com
Technical Support: tech@systembio.com

Ordering Information: orders@systembio.com

Licensing and Warranty Statement

Limited Use License

Use of the ExoGlow-NTA Fluorescent Labeling Kit (*i.e.*, the “Product”) is subject to the following terms and conditions. If the terms and conditions are not acceptable, return all components of the Product to System Biosciences (SBI) within 7 calendar days. Purchase and use of any part of the Product constitutes acceptance of the above terms.

The purchaser of the Product is granted a limited license to use the Product under the following terms and conditions:

- The Product shall be used by the purchaser for internal research purposes only. The Product is expressly not designed, intended, or warranted for use in humans or for therapeutic or diagnostic use.
- The Product may not be resold, modified for resale, or used to manufacture commercial products without prior written consent of SBI.
- This Product should be used in accordance with the NIH guidelines developed for recombinant DNA and genetic research.

SBI has pending patent applications related to the Product. For information concerning licenses for commercial use, contact SBI.

Purchase of the product does not grant any rights or license for use other than those explicitly listed in this Licensing and Warranty Statement. Use of the Product for any use other than described expressly herein may be covered by patents or subject to rights other than those mentioned. SBI disclaims any and all responsibility for injury or damage which may be caused by the failure of the buyer or any other person to use the Product in accordance with the terms and conditions outlined herein.

Limited Warranty

SBI warrants that the Product meets the specifications described in this manual. If it is proven to the satisfaction of SBI that the Product fails to meet these specifications, SBI will replace the Product or provide the purchaser with a refund. This limited warranty shall not extend to anyone other than the original purchaser of the Product. Notice of nonconforming products must be made to SBI within 30 days of receipt of the Product.

SBI’s liability is expressly limited to replacement of Product or a refund limited to the actual purchase price. SBI’s liability does not extend to any damages arising from use or improper use of the Product, or losses associated with the use of additional materials or reagents. This limited warranty is the sole and exclusive warranty. SBI does not provide any other warranties of any kind, expressed or implied, including the merchantability or fitness of the Product for a particular purpose.

SBI is committed to providing our customers with high-quality products. If you should have any questions or concerns about any SBI products, please contact us at (888) 266-5066.

This page intentionally left blank.

This page intentionally left blank.

This page intentionally left blank.

This page intentionally left blank.



System Biosciences (SBI)
2438 Embarcadero Way
Palo Alto, CA 94303

Phone: (650) 968-2200
Toll-Free: (888) 266-5066
Fax: (650) 968-2277

E-mail:
General Information: info@systembio.com
Technical Support: tech@systembio.com
Ordering Information: orders@systembio.com